Check Stab Instruments

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Functions of the Check Stab α 2012 Life

- *Quick and easy measurement of tartaric acid stability using electrochemical conductivity measurements at various conditions.*
- *Mini Contact: Determines KHT precipitation expressed in μ Siemens. This value indicates the relative stability of the wine sample being tested.*
- *Mini contact with forecast: Determines KHT precipitation expressed in μ Siemen and by using CheckFuture®, predict future precipitation of crystals.*
- *Saturation Point: Automatically, the analysis calculates the saturation temperature expressed in °C of the test sample. It has also the possibility to evaluate the data with the Total Polyphenols Index.*
- *Saturation Point of Calcium: This analysis calculates automatically the saturation temperature value of calcium.*

*Due to numerous requests from many of our customers, **CheckStab Instruments®**, leader in the development of instruments for the analysis of tartaric stability in wines, has developed a new generation instrument. The latest unit is the **Check Stab α 2012 Life**, a completely automatic analyzer capable of satisfying all types of wineries, small or large. The laboratory can connect from 1 to 16 instruments together, all controlled by one software CheckStab Net®, with the data stored on one single Data Base. The instrument has an automatic KHT dispensing system. The connection of the analyzer to the computer is done by means of LAN which makes data management much more simplified thanks to the standardization of the communication protocol.*

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View of the automatic KHT dispenser

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Some characteristics of the Check Stab α 2012 Life

The operation and management of data is done with a PC with a Windows operating system. The instruments can be connected to the PC by LAN, WLAN, USB or Serial Port Rs 232c. Possibility to connect to Bar Code reader to insert sample data. Possibility to connect more than one instrument together (max 16 units), with each unit doing an independent analysis. Automatic dispensing system for adding KHT without the intervention of the lab technician. Stand by function for energy savings.

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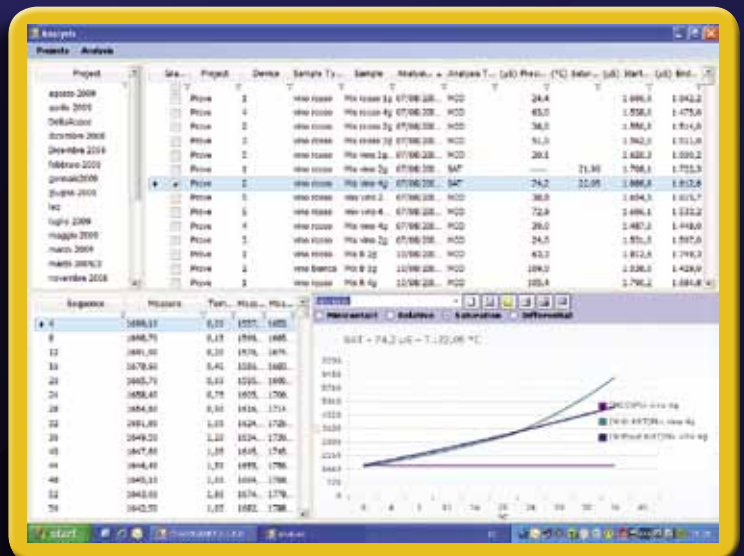
Acquiring and recording measurement data is becoming more and more important. CheckStab Instruments® has developed a new analytical software system to manage data. This permits the lab to study, store, and comment on an analysis with extreme simplicity. The entire system is managed by use of a personal computer.

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What Check Stab α 2012 Life can do

- *Contemporaneous management of connected instruments*
- *Possibility to do contemporaneously several analyses, with independent parameter research*
- *Visualize measurement data (Conductivity, Wine temperature, Bath temperature, pre-set temperature, etc.).*
- *Visualize user status, using incorporated graphic symbols (State of compressor, resistance, probe positions, etc.).*
- *Memorize the analysis on the data-base, with the possibility to transfer to a text file.*
- *Possibility to get from Check Stab analyses made, to visualize results in graphic form, to visualize results in text form, to save on a disk, export reports in various formats (PDF, Excel, Text) or print the graphics using a printer.*
- *Visualize and print the analysis in graphics or text format.*
- *Visualize the saturation analysis in a normal graphics mode or in a differential mode.*
- *Visualize the mini contact analysis in absolute mode or in relative mode.*
- *Visualize the analysis in single mode or multiple mode, up to a maximum of 10 analysis contemporaneously.*
- *Interface for LIMS*

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Data given in this brochure are indicative and can vary at the discretion of the manufacturer at any time and without obligation or notice.

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For local sales enquiries, please contact Vitis & Winemakers;

- Head Office: + 61 (03) 9487-1160
- NZ Office & Shop: +64 3 577-8778
- info@vitiswinemakers.com

Vitis & Winemakers www.vitiswinemakers.com